

Linearity Calibration

- Before linearity calibration, please adjust the level indicator firstly.
1. Please connect Port No. 5 and No. 6 of the RS232C interface. We will offer a special connector for you.
 2. Press the key "COU", it will show "LIN-200", then put on 200g weight, the balance will show "-----", when it show "LIN-0", take away the weight.
 3. Then the balance will show "LIN-100" automatically, please put 100g weight on the pan, it will show "-----", when it show "LIN-0", take away the weight.
 4. Now the balance will show "LIN-50", put on 50 g weight, then show "-----", when it show "LIN-0", take away the weight too.
 5. At last, the balance will show "LIN-150", put on 150g weight, then it show "-----", the balance will show "CAL-0-", then "CAL-200" flicker, please put 200g weight on the plate. When the display is stable, calibration is ok.

Please note, our all Magnetic Balance 0.1mg and 1mg have the same operation ways of linearity calibration. So just follow the instructions of the Balance LCD display. TKS.